



Full Range Pipette (1 - 1000  $\mu$ L)

---

**Quick Start Guide**

2-APR-2026

**DeNovix**<sup>®</sup>

## Active Pipetting Mode

See User Guide for pipetting mode instructions.


## Tip Status

## Action Wheel

Hold for 0.5 seconds to wake from Sleep Mode.

Press to aspirate, dispense or select.

Turn to scroll or adjust volume.

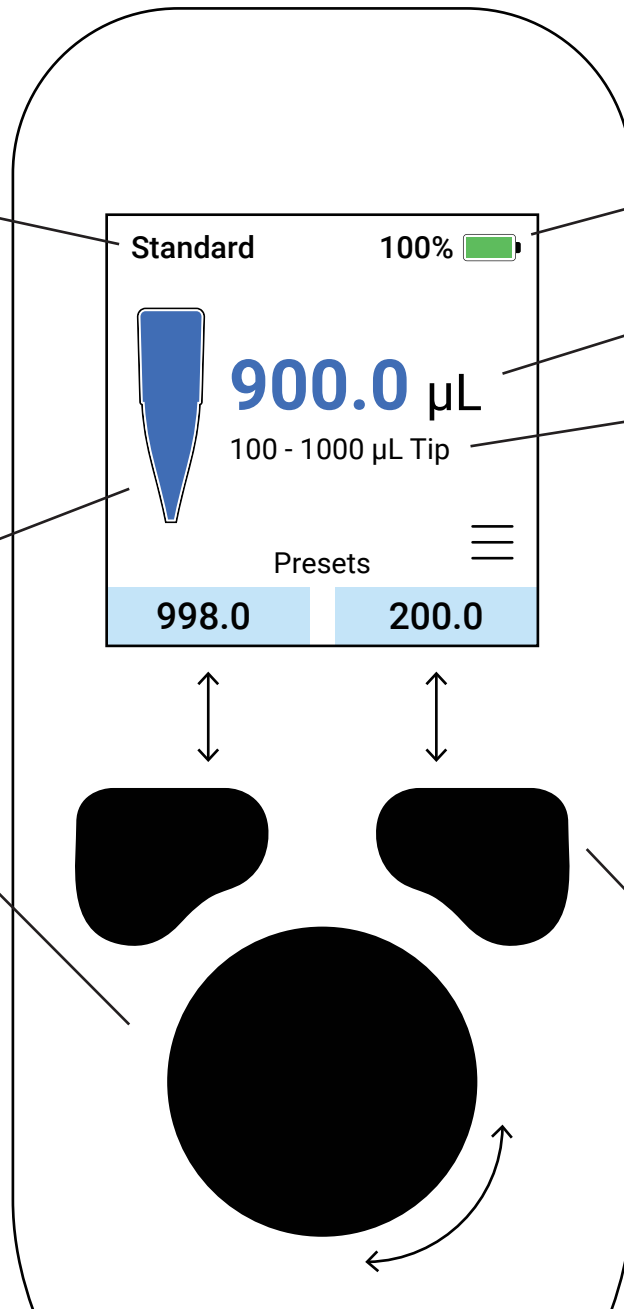

## Battery

## Volume

## Tip Mode

## Side Button

Access menu or return to previous screen.

Press and hold to put into Sleep Mode.

## Left and Right Buttons

Press to activate Presets.

Hold to edit value.

# Squid™ Full Range Pipette

Squid™ is a 1 - 1000 µL electronic pipette that uses two tip sizes for different volume ranges.

- Large Tip Mode: 100 - 1000 µL
- Small Tip Mode: 1 - 200 µL


## What's in the Box?

- Squid™ Full Range Pipette (1 - 1000 µL)
- Charging Cable and Charger
- Certification
- O-Ring Grease
- Extra O-Rings

## Getting Started

1. Remove Squid Pipette from packaging.
2. Hold Action Wheel for 0.5 seconds to wake from Sleep Mode.
3. Use the Side Button to select a pipetting mode or adjust settings.

*For detailed instructions, please refer to the User Guide (QR code on back).*

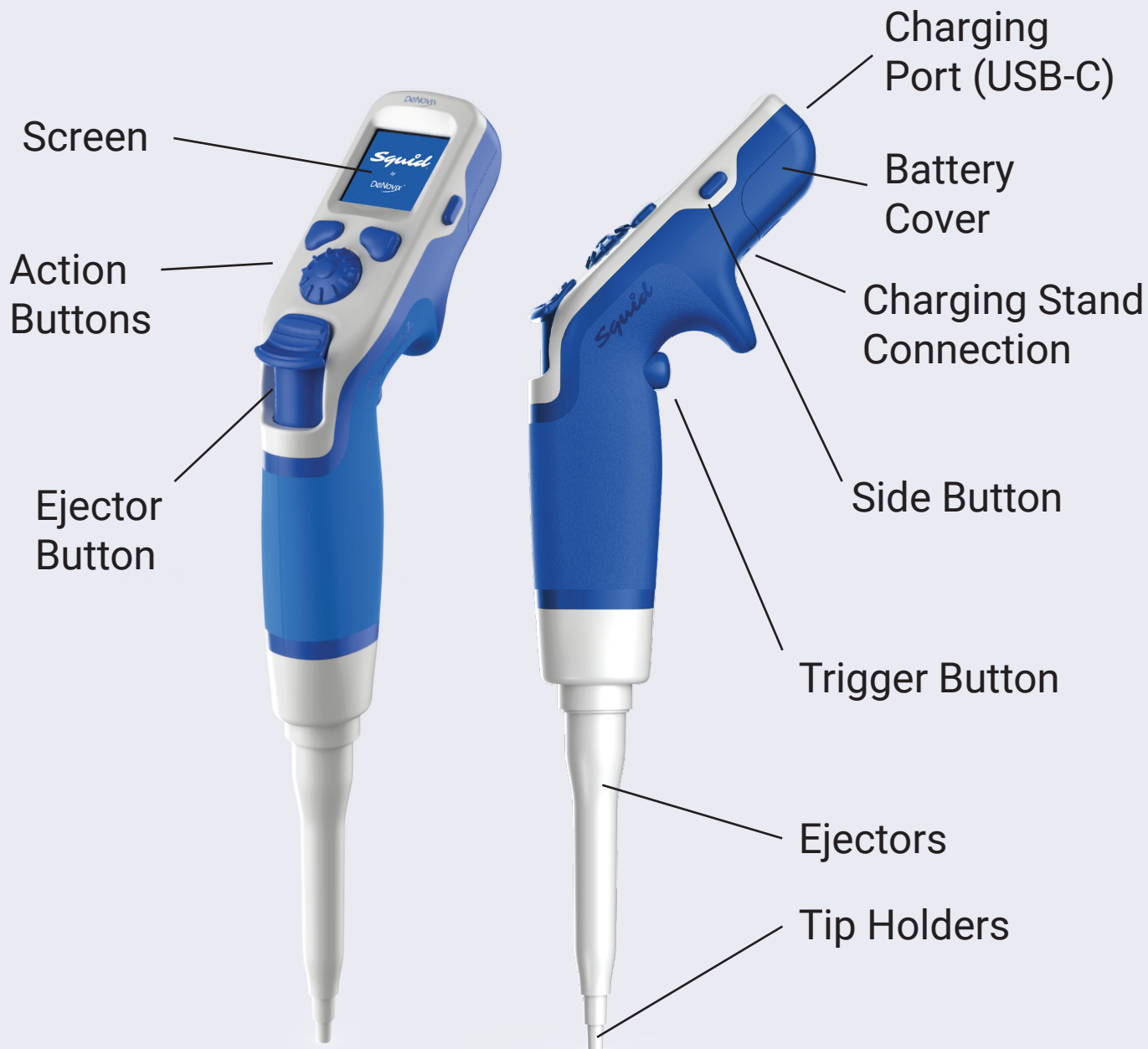

# Using Squid Tip Modes

## Large Tip Mode

100 - 1000  $\mu$ L



Press Ejector



Large Tip Ready

## Small Tip Mode

1 - 200  $\mu$ L




Press Trigger Button



Small Tip Ready

**Important:** After ejecting, press Trigger Button to return to Small Tip Mode.

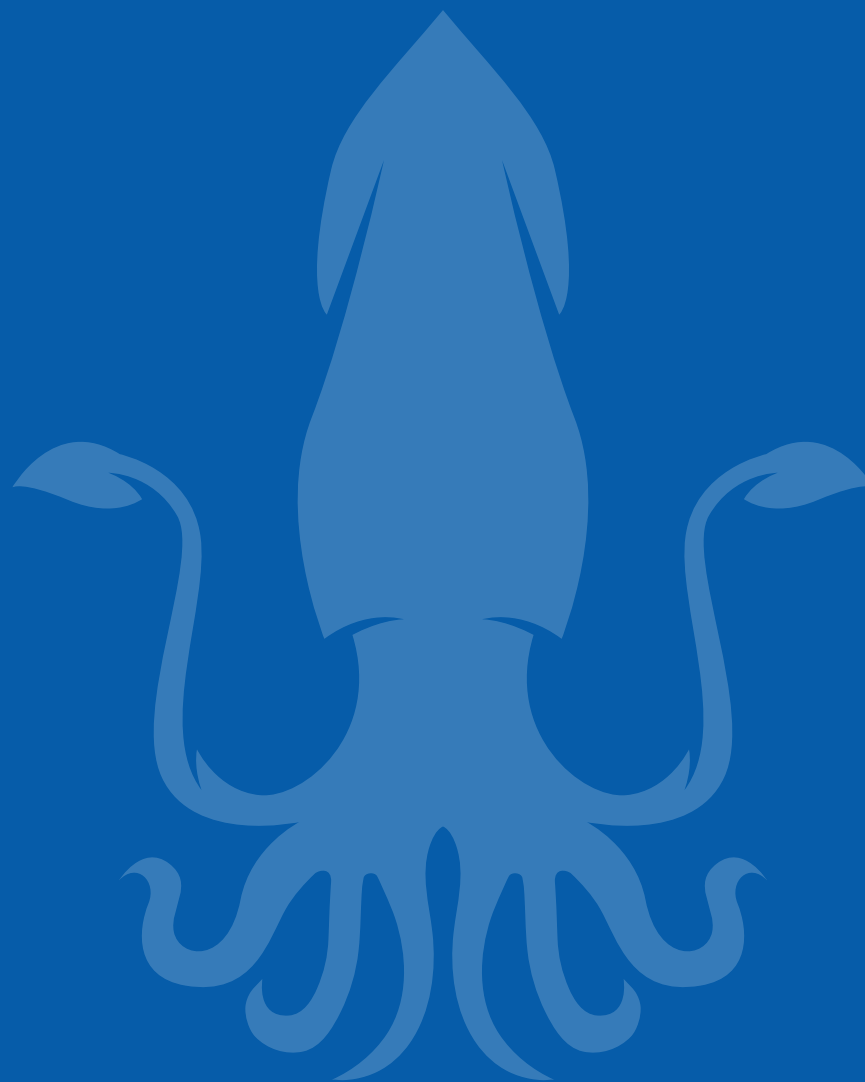

**Biozym**<sup>®</sup>

SCIENCE IS OUR BUSINESS

[www.biozym.com](http://www.biozym.com)

**Biozym Scientific GmbH**

Tel.: +49 (0) 5152 9020

Mail: [support@biozym.com](mailto:support@biozym.com)

**Biozym Biotech Trading GmbH**

Tel.: +43 (0) 1 334 0156 0

Mail: [support@biozym.com](mailto:support@biozym.com)

