

DuraPorter® Sample Carrier

Important Information

- The unit should be locked when carried with samples enclosed.
- The unit should be discarded if the locks or housing are broken or cracked. Follow your institution's policies for the disposal of biological hazardous materials.
- The unit will house most standard size 13 mm and 16 mm tube racks.
- The unit includes dividers for organizing bagged or other sample containers.
- The sample carrier is autoclavable (121° C for 15 minutes). Higher temperatures or autoclaving duration may damage the unit. The unit should be unlocked and open during autoclaving with the silicone seal left in place.
- The sample carrier should not be used with radioactive materials.