



PRACTICAL INNOVATION

Biozym
SCIENCE IS OUR BUSINESS

Ziath

DataPaq™ Single-Tube Scanner

Fast, small, "point-and-scan", with a 2-year guarantee

The Ziath DataPaq™ Single-Tube Scanner scans single tubes from Matrix, Nunc, ABgene, Micronic, Greiner, Wheaton and all other tube manufacturers. Simply hold the tube to the large scanner window and the code is instantly displayed.



Features

- Fast scanning: less than 1 second per tube
- Compatible with Matrix, ABgene, Nunc, REMP, Micronic, Greiner, Wheaton tubes and many more
- Instant read for any size or brand of tube
- Export data to any application: Excel, LIMS, etc.
- Operating systems: Windows XP, Vista, 7
- Simple installer; no dongles required
- Compact: 60 x 90 x 60mm (WxDxH)
- Durable construction - 2 year guarantee
- Free software upgrades for life
- Ziath's expertise and support come as standard



PRACTICAL INNOVATION

DataPaq™ Single Tube Scanner

Easy Set-Up and Easy To Use

The DataPaq™ Single Tube Scanner offers easy set-up, straight out of the box, with only a few minutes for configuration.

After a one-time automatic calibration, the scanner is ready to read Matrix, ABgene, Nunc, Micronic, Greiner, Wheaton and all other tubes on the market.


Our DataPaq™ software makes it simple to export to Excel, XML or text and scanned images can also be saved.

The large scan window makes it easy to present the tubes, even when wearing cryoprotective clothing and the weighted scanner base makes it a stable unit.

No independent power supply is required as the scanner connects to the computer by USB cable and also draws its power from this source

Cryoprotection™ Option

The occurrence of condensation whilst scanning tubes straight from cryogenic storage is well known. Ziath's cryoprotective coating eliminates the occurrence and ensures uninterrupted scanning. As Ziath's method for eliminating condensation uses no extra electrical components reliability is increased and costs are significantly decreased.

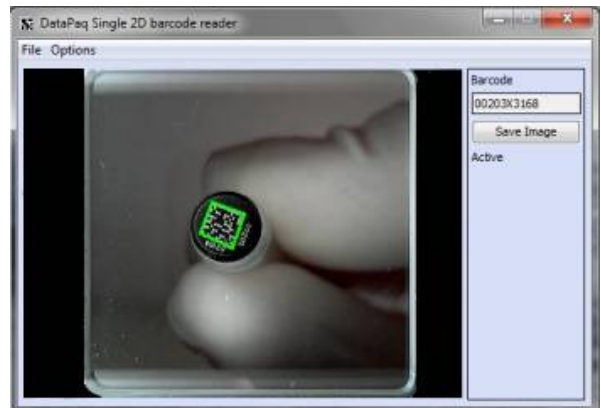

It's simple to export to Excel, XML or text using the DataPaq™ software



DataPaq Single Tube Scanner without Cryoprotection™. After repeated scanning, condensation starts to appear



DataPaq Single Tube Scanner with Cryoprotection™. No condensation has formed

Ordering Information

Code	Description	Quantity
601071	DataPaq™ Single Tube Scanner	1
601072	DataPaq™ Single Tube Scanner with Cryoprotection™	1

Biozym Scientific GmbH
Tel: +49 (0)5152-9020
Fax: +49 (0)5152-2070
support@biozym.com www.biozym.com

